碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301或800-8283301
订货e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

Ad-mCherry-GFP-LC3B

产品编号	产品名称	包装
C3011-1ml	Ad-mCherry-GFP-LC3B	1ml
C3011-10ml	Ad-mCherry-GFP-LC3B	10ml

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Species	Cell line	Origin	Transduction efficiency
Human	293	Embryonic Kidney	95%
	MCF-10A	Mammary epithelium	51%
	MDA-MB-231	Malignant breast cancer cell	72%
	HepG2	Hepatocarcinoma	70%
	T24	Transitional cell carcinoma	37%
	TCCSUP	Transitional cell carcinoma	53%
	PC-3	Prostate cancer metastases site	67%
	DU145	Prostate cancer metastases site	86%
	LNCaP	Prostate cancer metastases site	69%
	Jurkat	T cell leukemia	2%
	HRT	Rectal tumor	5%
Murine	MT1A2	Mammary tumor transformed with SV40 T antigen	92%
	NIH 3T3	Mouse fibroblast	12%
	WEHI	Fibrosarcoma	10%
	AML12	Liver	28%
	MS-K68	Spleen	48%
	EL4	T cell	2%
Bovine	FBRT HE1	Fetal retina transformed with Had5 E1	90%
	MDBK	Kidney	70%
	FBK-34	Kidney transformed with BAd3 E1	25%
	BT-5705	Turbinate	28%
	EBL	Embryonic lung	8%
Porcine	FPRT HE1-5	Fetal retina transformed with HA5 E1	91%
	PK-15	Kidney	76%
	PK-M12	Kidney cell transformed with SV40 T antigen	71%
	PT-K75	Turbinate	22%
	EPL	Embryonic lung	53%
	SBC-1765	Derived from buffy coat cells	73%

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Table 1. Summary of Recommended Biosafety Levels for Infectious Agents

BSL	Agents	Practices	Primary Barriers and Safety Equipment	Facilities (Secondary Barriers)
-----	--------	-----------	---------------------------------------	---------------------------------

1	Not known to consistently cause diseases in healthy adults	Standard microbiological practices	<ul style="list-style-type: none"> ■ No primary barriers required. ■ PPE: laboratory coats and gloves; eye, face protection, as needed 	Laboratory bench and sink required
2	<ul style="list-style-type: none"> ■ Agents associated with human disease ■ Routes of transmission include percutaneous injury, ingestion, mucous membrane exposure 	BSL-1 practice plus: <ul style="list-style-type: none"> ■ Limited access ■ Biohazard warning signs ■ “Sharps” precautions ■ Biosafety manual defining any needed waste decontamination or medical surveillance policies 	Primary barriers: <ul style="list-style-type: none"> ■ BSCs or other physical containment devices used for all manipulations of agents that cause splashes or aerosols of infectious materials ■ PPE: Laboratory coats, gloves, face and eye protection, as needed 	BSL-1 plus: <ul style="list-style-type: none"> ■ Autoclave available
3	Indigenous or exotic agents that may cause serious or potentially lethal disease through the inhalation route of exposure	BSL-2 practice plus: <ul style="list-style-type: none"> ■ Controlled access ■ Decontamination of all waste ■ Decontamination of laboratory clothing before laundering 	Primary barriers: <ul style="list-style-type: none"> ■ BSCs or other physical containment devices used for all open manipulations of agents ■ PPE: Protective laboratory clothing, gloves, face, eye and respiratory protection, as needed 	BSL-2 plus: <ul style="list-style-type: none"> ■ Physical separation from access corridors ■ Self-closing, double-door access ■ Exhausted air not recirculated ■ Negative airflow into laboratory ■ Entry through airlock or anteroom ■ Hand washing sink near laboratory exit
4	<ul style="list-style-type: none"> ■ Dangerous/exotic agents which pose high individual risk of aerosol-transmitted laboratory infections that are frequently fatal, for which there are no vaccines or treatments ■ Agents with a close or identical antigenic relationship to an agent requiring BSL-4 until data are available to redesignate the level ■ Related agents with unknown risk of transmission 	BSL-3 practices plus: <ul style="list-style-type: none"> ■ Clothing change before entering ■ Shower on exit ■ All material decontaminated on exit from facility 	Primary barriers: <ul style="list-style-type: none"> ■ All procedures conducted in Class III BSCs or Class I or II BSCs in combination with full-body, air-supplied, positive pressure suit 	BSL-3 plus: <ul style="list-style-type: none"> ■ Separate building or isolated zone ■ Dedicated supply and exhaust, vacuum, and decontamination systems ■ Other requirements outlined in the text

BSL, biosafety level; PPE, personal protective equipment.

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产品编号	产品名称	包装
C3006-1ml	Ad-GFP-LC3B	1ml
C3006-10ml	Ad-GFP-LC3B	10ml
C3011-1ml	Ad-mCherry-GFP-LC3B	1ml
C3011-10ml	Ad-mCherry-GFP-LC3B	10ml
C3012-1ml	AdPlus-mCherry-GFP-LC3B	1ml
C3012-10ml	AdPlus-mCherry-GFP-LC3B	10ml

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1. Yu P, Zhang C, Gao CY, Ma T, Zhang H, Zhou MM, Yang YW, Yang L, Kong LY. Anti-proliferation of triple-negative breast cancer cells with physagulide P: ROS/JNK signaling pathway induces apoptosis and autophagic cell death. *ONCOTARGET* . 2017 Jul 17;8(38):64032-64049.
2. Peng Y, Qiu L, Xu D, Zhang L, Yu H, Ding Y, Deng L, Lin J. M4IDP, a zoledronic acid derivative, induces G1 arrest, apoptosis and autophagy in HCT116 colon carcinoma cells via blocking PI3K/Akt/mTOR pathway. *Life Sci* . 2017 Sep 15;185:63-72.
3. Ma Z, Wang C, Liu C, Yan DY, Deng Y, Liu W, Yang TY, Xu ZF, Xu B. The role S-nitrosylation in manganese-induced autophagy dysregulation in SH-SY5Y cells. *Environ Toxicol* . 2017 Dec;32(12):2428-2439.
4. He X, Yuan W, Li Z, Hou Y, Liu F, Feng J. 6-Hydroxydopamine induces autophagic flux dysfunction by impairing transcription factor EB activation and lysosomal function in dopaminergic neurons and SH-SY5Y cells. *Toxicol Lett* . 2018 Feb;283:58-68.
5. He X, Yuan W, Li Z, Hou Y, Liu F, Feng J. 6-Hydroxydopamine induces autophagic flux dysfunction by impairing transcription factor EB activation and lysosomal function in dopaminergic neurons and SH-SY5Y cells. *Toxicol Lett* . 2018 Feb;283:58-68.
6. Xiao Y, Zhou Y, Lu Y, Zhou K, Cai W. PHB2 interacts with LC3 and SQSTM1 is required for bile acids-induced mitophagy in cholestatic liver. *Cell Death Dis*. 2018 Feb 7;9(2):160.

7. Wang J,Hussain T,Yue R,Liao Y,Li Q,Yao J,Song Y,Sun X,Wang N,Xu L,Sreevatsan S,Zhao D,Zhou X.MicroRNA-199a Inhibits Cellular Autophagy and Downregulates IFN- β Expression by Targeting TBK1 in Mycobacterium bovis Infected Cells.Front Cell Infect Microbiol . 2018 Jul 10;8:238.
8. Zhang Z,Wei Y,Liu B,Wu Y,Wang H,Xie X,Feng Z,Shao G,Xiong Q.Hsp90/Sec22b promotes unconventional secretion of mature-IL-1 β through an autophagosomal carrier in porcine alveolar macrophages during Mycoplasma hyopneumoniae infection.Mol Immunol . 2018 Sep;101:130-139.
9. Liu M,Zhao G,Zhang D,An W,Lai H,Li X,Cao S,Lin X.Active fraction of clove induces apoptosis via PI3K/Akt/mTOR-mediated autophagy in human colorectal cancer HCT-116 cells.Int J Oncol . 2018 Sep;53(3):1363-1373.
10. Yin J,Ni B,Yang YD,Tang ZW,Gao ZQ,Feng L,Liao WG,Gao YQ.Elevation of autophagy rescues spermatogenesis by inhibiting apoptosis of mouse spermatocytes.Reproduction . 2018 Sep 27. pii: REP-18-0243.
11. Yang Y,Gao J,Zhang Y,Xu W,Hao Y,Xu Z,Tao L.Natural pyrethrins induce autophagy of HepG2 cells through the activation of AMPK/mTOR pathway.Environ Pollut . 2018 Oct;241:1091-1097.
12. Wei R,Mao L,Xu P,Zheng X,Hackman RM,Mackenzie GG,Wang Y.Suppressing glucose metabolism with epigallocatechin-3-gallate (EGCG) reduces breast cancer cell growth in preclinical models.Food Funct . 2018 Nov 14;9(11):5682-5696.
13. Zhuo FF,Zhang C,Zhang H,Xia Y,Xue GM,Yang L,Kong LY.Chrysanthemulide A induces apoptosis through DR5 upregulation via JNK-mediated autophagosome accumulation in human osteosarcoma cells.J Cell Physiol . 2018 Dec 17.
14. Liu B,Ding L,Zhang L,Wang S,Wang Y,Wang B,Li L.Baicalein Induces Autophagy and Apoptosis through AMPK Pathway in Human Glioma Cells.Am J Chin Med. 2019;47(6):1405-1418.
15. Ma Z,Li F,Chen L,Gu T,Zhang Q,Qu Y,Xu M,Cai X,Lu L.Autophagy promotes hepatic differentiation of hepatic progenitor cells by regulating the Wnt/ β -catenin signaling pathway.J Mol Histol. 2019 Feb;50(1):75-90.
16. Yan D,Ma Z,Liu C,Wang C,Deng Y,Liu W,Xu B.Corynoxine B ameliorates HMGB1-dependent autophagy dysfunction during manganese exposure in SH-SY5Y human neuroblastoma cells.Food Chem Toxicol. 2019 Feb;124:336-348.
17. Yang X,Gao J,Guo J,Zhao Z,Zhang SL,He Y.Anti-lung cancer activity and inhibitory mechanisms of a novel Calothrixin A derivative.Life Sci. 2019 Feb 15;219:20-30.
18. Chu Q,Zhang S,Chen M,Han W,Jia R,Chen W,Zheng X.Cherry Anthocyanins Regulate NAFLD by Promoting Autophagy Pathway.Oxid Med Cell Longev. 2019 Feb 25;2019:4825949.
19. Luo Z,Xu X,Sho T,Zhang J,Xu W,Yao J,Xu J.ROS-induced autophagy regulates porcine trophectoderm cell apoptosis, proliferation, and differentiation.AM J PHYSIOL-CELL PH. 2019 Feb 1;316(2):C198-C209.
20. Ren Z,Xiao W,Zeng Y,Liu MH,Li GH,Tang ZH,Qu SL,Hao YM,Yuan HQ,Jiang ZS.Fibroblast growth factor-21 alleviates hypoxia/reoxygenation injury in H9c2 cardiomyocytes by promoting autophagic flux.Int J Mol Med. 2019 Mar;43(3):1321-1330.
21. Wang X,Xiao D,Ma C,Zhang L,Duan Q,Zheng X,Mao M,Zhu D,Li Q.The effect of honokiol on pulmonary artery endothelium cell autophagy mediated by cyclophilin A in hypoxic pulmonary arterial hypertension.J Pharmacol Sci. 2019 Mar;139(3):158-165.
22. Zeng H,Zhang W,Gong Y,Xie C.Radiotherapy activates autophagy to increase CD8+ T cell infiltration by modulating major histocompatibility complex class-I expression in non-small cell lung cancer.J Int Med Res. 2019 Aug;47(8):3818-3830.
23. Gao S,Li S,Li Q,Zhang F,Sun M,Wan Z,Wang S.Protective effects of salvianolic acid B against hydrogen peroxide-induced apoptosis of human umbilical vein endothelial cells and underlying mechanisms.Int J Mol Med. 2019 Aug;44(2):457-468.
24. Chu P,He L,Yang C,Zeng W,Huang R,Liao L,Li Y,Zhu Z,Wang Y.Grass carp ATG5 and ATG12 promote autophagy but down-regulate the transcriptional expression levels of IFN- γ signaling pathway.Fish Shellfish Immunol. 2019 Sep;92:600-611.
25. Huang Y,Li Y,Qu Y,Zheng Y,Ouyang M,Zhang Y,Lai W,Xu Q.UVA-induced photodamage inhibits autophagic degradation by impairing lysosomal function in dermal fibroblasts.Biochem Biophys Res Commun. 2019 Oct 22;518(4):611-618.
26. Li R,Du J,Yao Y,Yao G,Wang X.Adiponectin inhibits high glucose-induced angiogenesis via inhibiting autophagy in RF/6A cells.J Cell Physiol. 2019 Nov;234(11):20566-20576.
27. Hong X,Min SN,Zhang YY,Lin YT,Wang F,Huang Y,Yu GY,Wu LL,Yang HY.TNF- α Suppresses Autophagic Flux in Acinar Cells in IgG4-Related Sialadenitis.J Dent Res. 2019 Nov;98(12):1386-1396.
28. He J,Shi XY,Li ZM,Pan XH,Li ZL,Chen Y,Yan SJ,Xiao L.Proton pump inhibitors can reverse the YAP mediated paclitaxel resistance in epithelial ovarian cancer.BMC Mol Cell Biol. 2019 Nov 12;20(1):49.
29. Zhang Y,Guo W,Chen H,Gao J,Xu Z,Tao L,Li Z,Xu W.Spinetoram confers its cytotoxic effects by inducing AMPK/mTOR-mediated autophagy and oxidative DNA damage.Ecotoxicol Environ Saf. 2019 Nov 15;183:109480.
30. Yan Z,Su G,Gao W,He J,Shen Y,Zeng Y,Liu X.Fluid shear stress induces cell migration and invasion via activating autophagy in HepG2 cells.Cell Adh Migr. 2019 Dec;13(1):152-163.
31. Zhuo Ma,Can Wang,Chang Liu,Dong-Ying Yan,Xuan Tan,Kuan Liu,Meng-Jiao Jing,Yu Deng,Wei Liu,Bin Xu.Manganese induces autophagy dysregulation: The role of S-nitrosylation in regulating autophagy related proteins in vivo and in vitro.Sci Total Environ. 2020 Jan 1;698:134294.;doi: 10.1016/j.scitotenv.2019.134294
32. Chang Liu,Dong-Ying Yan,Can Wang,Zhuo Ma,Yu Deng,Wei Liu,Bin Xu.Manganese activates autophagy to alleviate endoplasmic reticulum stress-induced apoptosis via PERK pathway J Cell Mol Med. 2020 Jan;24(1):328-341.;doi: 10.1111/jcmm.14732
33. Su Li,Zhilei Bo,Ying Jiang,Xianmin Song,Chun Wang,Yin Tong.Homoharringtonine promotes BCR-ABL degradation through the p62-mediated autophagy pathway Oncol Rep. 2020 Jan;43(1):113-120.;doi: 10.3892/or.2019.7412
34. Chang Liu,Dong-Ying Yan,Can Wang,Zhuo Ma,Yu Deng,Wei Liu,Bin Xu.IRE1 signaling pathway mediates protective autophagic response against manganese-induced neuronal apoptosis in vivo and in vitro Sci Total Environ. 2020 Apr 10;712:136480.;doi: 10.1016/j.scitotenv.2019.136480
35. Na Geng,Xiaozhou Wang,Xiaohui Yu,Run Wang,Yiran Zhu,Meihua Zhang,Jianzhu Liu,Yongxia Liu.Staphylococcus aureus Avoids Autophagy Clearance of Bovine Mammary Epithelial Cells by Impairing Lysosomal Function Front Immunol. 2020 May 5;11:746.;doi: 10.3389/fimmu.2020.00746
36. Ying Wang,Xin Zhang,Wenjia Chen,Lei Gao,Jihe Li,Tao Song,Jinyu Chi,Xiaohui Zhang,Zhiyu Shi,Yanghong Dong,Xinhua Yin,Yue Liu.Cortistatin ameliorates Ang II-induced proliferation of vascular smooth muscle cells by inhibiting autophagy through SSTR3 and SSTR5 Life Sci. 2020 Jul 15;253:117726.;doi: 10.1016/j.lfs.2020.117726
37. Chunli Xu,Qingke Bai,Chen Wang,Qiuyu Meng,Yuming Gu,Qiwei Wang,Wenjia Xu,Ying Han,Yong Qin,Song Jia,Junfang Zhang,Jie Xu,Jiao Li,Miao Chen,Feng Wang.miR-433 Inhibits Neuronal Growth and Promotes Autophagy in Mouse Hippocampal HT-22 Cell Line Front Pharmacol. 2020 Nov 23;11:536913.;doi: 10.3389/fphar.2020.536913
38. Yong Zhao,Minling Zhu,Wenjia Chen,Qing Chang,Li Shen,Runan Yan,Liuping Yang,Jiashu Li,Rui Xu,Yu Fu,Yue Liu.TFPI α alleviated vascular endothelial cell injury by inhibiting autophagy and the class III PI3K/Beclin-1 pathway Thromb Res. 2020 Nov;195:151-157.;doi: 10.1016/j.thromres.2020.07.017. (IF 2.869)

39. Lixin Zhang,Cui Ma,Xiaoying Wang,June Bai,Siyu He,Junting Zhang,Wei Xin,Yiying Li,Yuan Jiang,Jianshuang Li,Daling Zhu.MicroRNA-874-5p regulates autophagy and proliferation in pulmonary artery smooth muscle cells by targeting Sirtuin 3 Eur J Pharmacol. 2020 Dec 5;888:173485.;doi: 10.1016/j.ejphar.2020.173485)
40. Li Zhang,Shixia Zhou,Tiejun Zhou,Kaifeng Yuan,Xiaoming Li,Junling Tang.S100A8 promotes chemoresistance via augmenting autophagy in B-cell lymphoma cells Oncol Rep. 2021 Jan;45(1):151-158.;doi: 10.3892/or.2020.7841
41. Li-Juan Gao,Ping Li,Tengyun Ma,Zhan-Qiong Zhong,Shi-Jun Xu.Ligustilide alleviates neurotoxicity in SH-SY5Y cells induced by A β 25-35 via regulating endoplasmic reticulum stress and autophagy Phytother Res. 2021 Mar;35(3):1572-1584.;doi: 10.1002/ptr.6925

Version 2021.09.01